


Addendum to MiniScan EZ User's Manual

Description

This CMR provides a 11-0050-30 Black Glass, 4-in x 4-in x 6-mm, uncalibrated, and a special 100-mm x 100-mm (4-in x 4-in) holder in which the regular MiniScan EZ 4500L Calibrated White Tile standards is placed for instrument standardization.

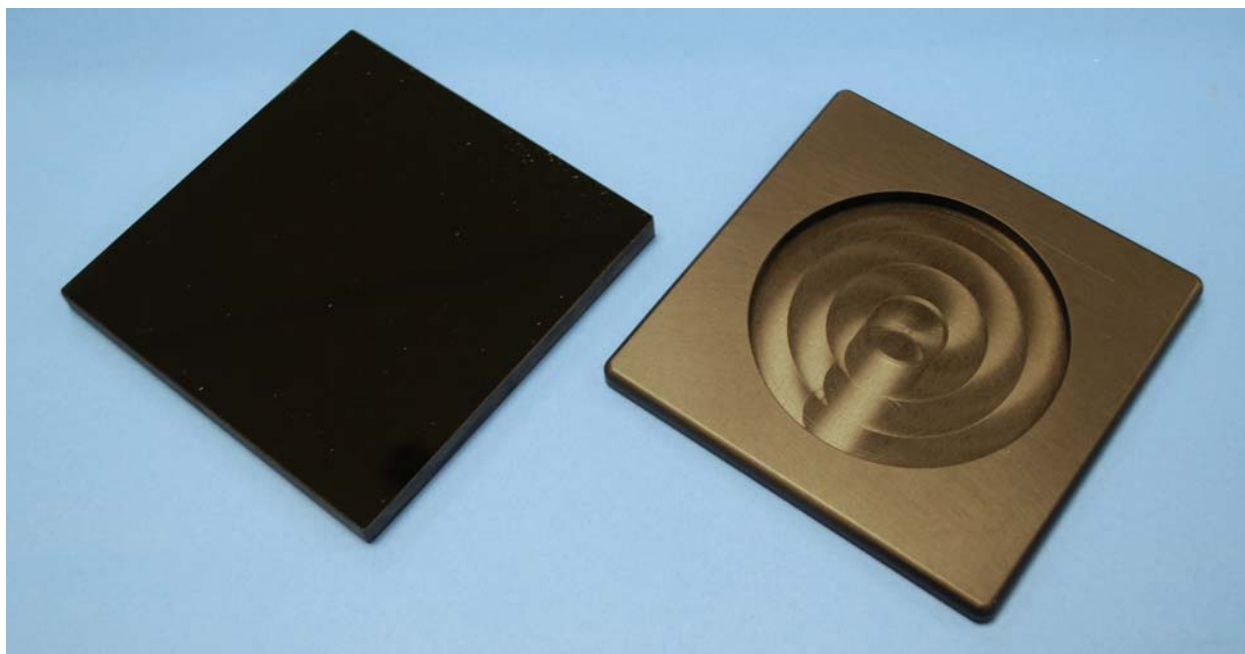

All MiniScan EZ replacement orders for Owens Corning should include the following items:

1. MSEZ 45/0 LAV instrument with standard options
2. D02-1014-367 – Nose Cone with Lower Glass Assembly to minimize dust in the instrument. This can be user-installed but should have been installed at the factory with a new instrument.
3. CMR3059 MSEZ 45/0 LAV 4-in x 4-in x 6-mm White Tile Holder + Black Glass to facilitate use with Owens Corning Standards Box

Installation

The main function of this CMR is to allow the use of the Owens Corning to use their Standards Box with 3 - 4-in x 4-in cutouts to hold the MSEZ Calibrated White Tile, a 4-in x 4-in x 6-mm Black Glass and a 4-in x 4-in OC Pink Tile (not part of this CMR).



All OC plants and affiliates are encouraged to communicate directly with Mike Stair to obtain the Owens Corning Standards Box (shown below) and A11-1008-596 Tile, 100-mm x 100-mm (4-in x 4-in), Pink Panther Tile. Contact info is:

Mr. Mike Stair
Owens Corning Corporation
Product Test Lab
2790 Columbus Road
Granville, OH USA 43023
T: (740) 321-7053
E: Mike.JMS.Stair@owenscorning.com

Operation

In the Owens Corning Standards Box, place the Black Glass in the far left slot and the MSEZ Calibrated White Tile from the standards cylinder placed in the CMR3059 Holder in the center. The OC Pink Panther Tile should be on the right.

Standardize normally using the Black Glass and Calibrated White Tile.

Measure the Pink Panther Tile as an OQ (Operation Qualification) step following the Owens Corning SOP.