


Addendum to MiniScan Diffuse User's Manual

Description

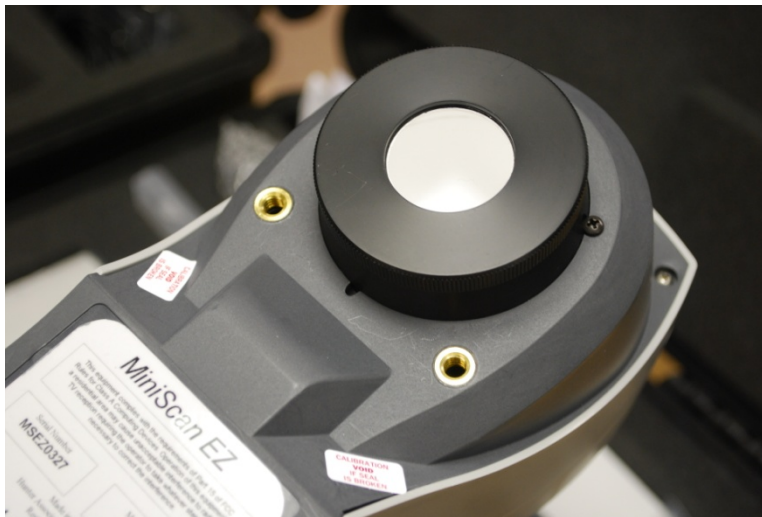

This CMR adds a field-replaceable glass port assembly to the MSEZ-4000L to protect the unit from dust. Also included are 5 extra pieces of replacement glass (A04-1001-920, Cover Glass) that can be replaced if the one currently in use becomes significantly scratched or abraded. For replacement, just unscrew the plastic faceplate and the glass port will pop out.

Installation

The instrument will come from the HunterLab factory with a special base that allows for a screw-on end plate with a round cover glass. The unit is ready for immediate operation.


If replacement glass is needed, contact HunterLab and order CMR3083-P3 Glass. To replace, unscrew the end cap and remove any parts of the old cover glass. Drop the new cover glass into the indentation of the end cap. Screw the end cap onto the base and the instrument is ready.

Operation

For taking normal color measurements, just standardize on the provided Light Trap and calibrated White Tile standards, then proceed to take readings.

Please note that instrument Diagnostic Tests must be run with an open port.

To perform White Tile short term repeatability and Green Tile long term diagnostic tests, the glass must first be removed and the end cap replaced such that there is an open port to match the conditions of assignment for these tests.

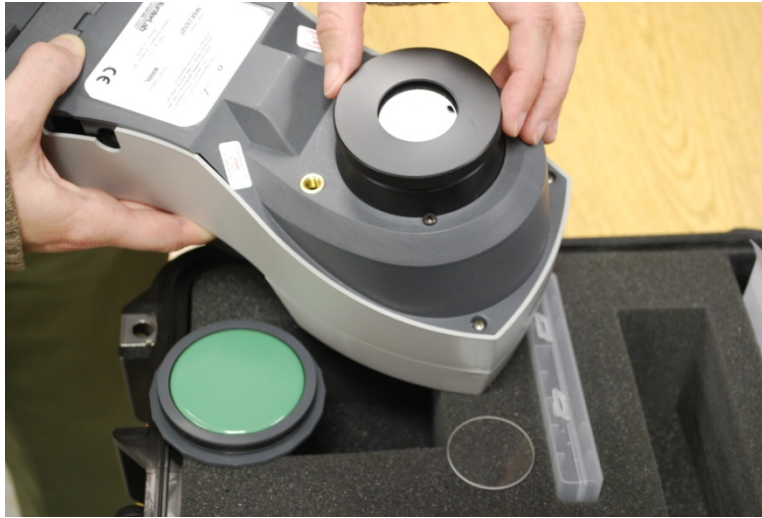

When finished with the diagnostic tests, replace the glass and standardize the instrument prior to making sample measurements.