

Values you Measured from Sample

Measured on Existing Instrument

Measured on Aeros

L or L*	a or a*	b or b*


Select **Numeric** Standard Type

Name Your Standard

Enter Values

Existing Standard or Target Value

L or L*	a or a*	b or b*

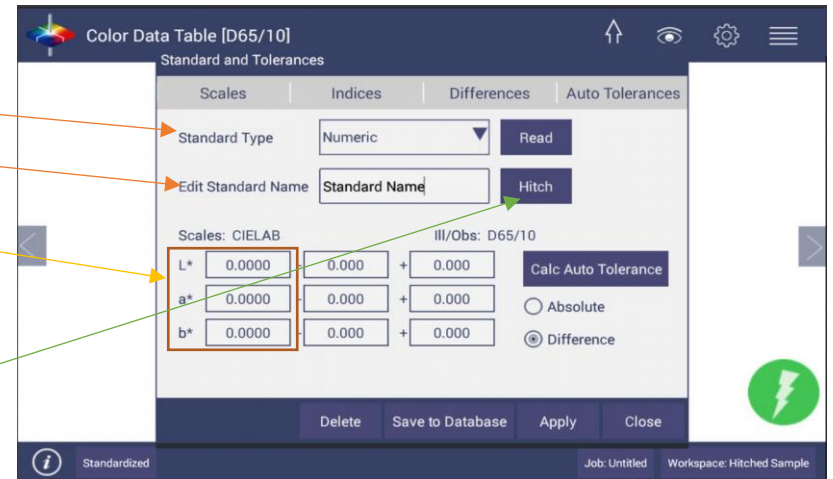


Press the Hitch Button

Enter these Calculated values into Target

0	0	0
---	---	---

Click Apply Hitch and then click OK

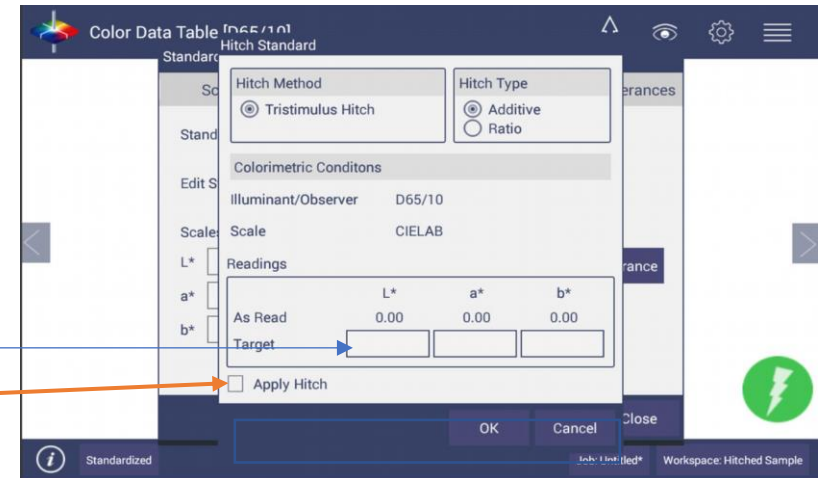

To make sure it is working correctly, re-measure the sample on Aeros. The values should match the Existing Instrument now, and the Standard shown in the Color Data Table should be your Existing Standard